

**Product:** **Chicken Embryo Extract**  
**Cat. No. 100-163P**  
Powder

---

#### Product Description

Gemini's chicken embryo extract is prepared from 9-12 day old eggs. It is processed as a 50% extract in phosphate-buffered saline that is supplemented with antibiotic:antimycotic. The liquid extract is sterile filtered to 0.2  $\mu\text{m}$  and lyophilized in 10mL aliquots. The sterility of this lyophilized powder cannot be guaranteed and should be sterile filtered once more after reconstitution.

#### Storage

Recommended storage is  $-20^{\circ}\text{C}$  or below. Lyophilized product may be stored at  $+4^{\circ}\text{C}$ .

#### Directions for Use

Reconstitute lyophilized product to 10ml using sterile water, agitate to mix.

It is advised that lyophilised product is filtered through 0.2 $\mu\text{m}$  filter after reconstitution to maintain sterility.

For *in vitro* diagnostic laboratory use or further manufacturing only. Not for human therapeutic use.